



# Automated Solid Phase Extraction of Endocrine Disruptors in Water

## Introduction

Endocrine disruptors need to be checked in drinking water at trace levels because of the effect they can have on hormones if allowed to reach high concentrations. SPE provides a quick and simple means of concentrating the components using controlled flow rates and volumes that will consistently produce good recoveries.

## Instrumentation Used for Sample Preparation



AUTOTRACE FOR AUTOMATED SPE AND THE TURBOVAP LV FOR EVAPORATION

## Sample Preparation

A 500ml sample of water is taken for the analysis of the following compounds.

- The solvent MTBE (Methylterbutylethylene ) is used as an Eluent.
- **Screen for Octylphenol, Nonylphenol, Octachlorostyrene, Butylbenzylphthalate, Diazinon, BHA & BHT.**
- The Eluent can be dried using an in line anhydrous sodium sulphate cartridge.

## AutoTrace Method

All solvent lines are purged & primed with solvent first. A maximum of 5 solvents can be used to run a wide range of methods. Independent lines separate the aqueous and organic waste solvents. The instrument sample rack has 6 positions that can take volumes from 10 ml to 2000 ml and 6 sample collection positions for the Eluent. This method offers an automated rugged and reproducible solution for cleaning up the samples to concentrate & remove interferences.

## AutoTrace Method

No.	Method:	Estimated time
		<b>1hr : 44 min</b>
1	Process 6 Samples using the following method steps:	
2	Condition the Column with 5.0 ml of Dichloromechane into solvent waste	
3	Condition Column with 5.0 ml of MTBE into solvent waste	
4	Condition Column with 5.0 ml of Methanol into solvent waste	
5	Condition Column with 5.0 ml of DI Water into aqueous waste	
6	Load 500.0 ml of sample onto Column	
7	Rinse Column with 5.0 ml of DI Water into aqueous waste	
8	Wash Syringe with 5.0 ml of MEOH/MTBE 10:90	
9	Dry Column with gas for 15.0 minutes	
10	Collect 5.0 ml Fraction into sample tube using MEOH/MTBE 10:90	
11	Collect 5.0 ml Faction into sample tube using Dichloromethane	
12	Wash Syringe with 5.0 ml of Methanol	
13	End	

### Parameters:

#### Flow Rates

Cond Flow:	15.0 ml/min
Load Flow:	10.0 ml/min
Rinse Flow:	20.0 ml/min
Elute Flow:	5.0 ml/min
Cond Air Push:	15.0 ml/min
Rinse Air Push:	20.0 ml/min
Elute Air Push:	5.0 ml/min

#### SPE Parameters

Push Delay:	5 sec
Air Factor:	1.0
Autowash Vol.:	1.00 ml

#### Workstation Parameters

Max. Elution Vol.:	12.0 ml
Exhaust Fan On:	Yes
Beeper On:	Yes



**Corporate Headquarters**  
68 Elm Street  
Hopkinton, MA 01748-1668

Tel: 508 436 9500  
Fax: 508 435 3439  
Email: [cust.support@caliperLS.com](mailto:cust.support@caliperLS.com)

[www.caliperLS.com](http://www.caliperLS.com)

©2008 Caliper Life Sciences. All rights reserved. Caliper, the Caliper logo, AutoTrace, RapidTrace, and TurboVap are trademarks or registered trademarks of Caliper Life Sciences. All other trademarks contained herein are the properties of their respective owners.

AT-AP-801 Mar 08